

CAY

For 1 liter of CAY:

1.5 g of casamino acids.

2.5 g of yeast extract.

20 ml of glycerol.

1 g of NaCl.

1 g of K_2HPO_4 .

0.1 g of $MgSO_4 \cdot 7H_2O$.

0.005 g of $CaCO_3$.

Complete to 1 liter with water.

For CAY-agar:

Add 15 g of agar per liter.

For CAY-top agar:

Add 7.5 g of agar per liter.

Autoclave.

Last modifications: 2011-01-24