

TYA

For 1 liter of TYA:

6 g of Tryptone.

2 g of Yeast extract.

3 g of Ammonium acetate.

0.2 g of $\text{MgSO}_4 \cdot 7\text{H}_2\text{O}$.

0.01 g of $\text{FeSO}_4 \cdot 7\text{H}_2\text{O}$.

Complete to 1 liter with water.

Autoclave.

Add 1% of sterile glucose and 3.7 mM of KH_2PO_4 .

For TYA-agar:

Add 15 g of agar per liter.

For TYA-top agar:

Add 7.5 g of agar per liter.

Autoclave.

Last modifications: 2011-02-25