

Media recipe

Czapeks-Dox + 0,2% YE

Détails

For 1 liter of Czapeks-Dox:

30 g of sucrose.

3 g of sodium nitrate.

1 g of dipotassium phosphate.

0.5 g of magnesium sulfate.

0.5 g of potassium chloride.

0.01 g of ferrous sulfate.

2 g of yeast extract.

Complete to 1 liter with water.

For Czapeks-Dox-agar:

Add 15 g of agar per liter.

For Czapeks-Dox-top agar:

Add 7.5 g of agar per liter.

Autoclave.

Last modifications: 2011-01-20