

Media recipe

Halobacterium broth (Halo)

Détails

For 1 liter of Halo:

250 g of NaCl.

20 g of $\text{MgSO}_4 \cdot 7\text{H}_2\text{O}$.

3 g of trisodium citrate $2\text{H}_2\text{O}$.

2 g of KCl.

5 g of yeast extract.

5 g of casamino acids.

2 g of L-arginine.

50 ml of Tris-HCl, 1M, pH7.4.

Complete to 1 liter with water.

For Halo-agar:

Add 15 g of agar per liter of 20% diluted broth.

Autoclave.

Last modifications: 2011-02-03