

Media recipe

Halobacterium cutirubrum medium (3.5S)

Détails

For 1 liter of 3.5S:

203 g of NaCl.

18 g of MgSO₄.

3.7 g of KCl.

0.11 g of CaCl₂.

1.26 mg of MnCl₂.

6 g of Yeast extract.

4 g of Casamino acids.

Complete to 1 liter with water.

For 3.5S-agar:

Add 15 g of agar per liter.

For 3.5S-top agar:

Add 7.5 g of agar per liter.

Autoclave.

Last modifications: 2011-02-19