

Media recipe

Alcalophilus medium (pH10)

Détails

For 1 liter of broth:

Part A:

900 ml of water
10 g of KH_2PO_4
5 g of peptone
5 g of yeast extract
1 g of glucose
0.2 g of $\text{MgSO}_4 \cdot 7\text{H}_2\text{O}$

Part B:

100 ml of water
20 g of Na_2CO_3

For agar:

Add 15 g of agar per liter to the part A

For top agar:

Add 7.5 g of agar per liter to the part A

Autoclave.

Mix the part A and part B.

Slepecky, R. A., and H. E. Hemphill. 1992. The Genus *Bacillus* - nonmedical, p. 1663-1696. In A.

B. e. al. (ed.), *The Prokaryotes*, 2nd ed. Springer-Verlag, New York Inc., Secaucus, New Jersey.

Dernières modifications: 16 janvier 2011