

Media recipe

Serum dextrose agar (SDA)

Détails

For 1 liter of SDA:

33 g of Difco Tryptose blood agar base	or	10 g of Tryptose
		3 g of Beef extract
		5 g of NaCl
		15 g of Agar

Complete to 950 ml with water.

Autoclave.

Add 50 ml of sterile horse serum.

Add 50 ml g of sterile glucose 20%.

Reference:

Corbel MJ, Thomas EL. 1980.

The Brucella-phages: their properties, characterization and applications. Booklet 2266, Ministry of Agriculture, Fisheries and Food, Pinner, Middlesex, UK, 106 p

Last modifications: 2011-02-24